amson Instruments Specification

Master Viscometers and Viscosity Oil Standards

ASTM D2162



Φ	Ultra-stable ± 0.005°C
\wedge	Intograted easiling

Integrated cooling Stainless steel bath

Heated window

RS232 communication

0.001°C readout

Item	Unit	TV16000	
P/N 230V/50Hz		00T0490	
P/N 115V/60Hz		00T0492	
P/N 230V/60Hz		00T0495	
Power	[kW]	3	
Range	[°C]	10 60	
Reading		°C or °F	
Setting	[°C]	0.001	
Stability ±	[°C]	< 0.005	
Uniformity ±	[°C]	< 0.005	
Heating	[kW]	Control: 1 + Boost: 1.5	
Heating total	[kW]	2.5	
Bath volume	[L]	160	
Bath opening	[mm]	250 x 365	
Bath depth	[mm]	965	
Length	[mm]	575	
Width	[mm]	575	
Height	[mm]	1700	
Weight	[kg]	135	
Materials	Used ins	Used inside bath: stainless steel 304, brass	
CE	Conforms to CE regulation		

General

This thermostatic bath is specially designed for tests that require ultra-precise temperature control, or processes that need to be followed visually. The TV16000 conforms to ASTM D2162: primary basic calibration of master viscometers and viscosity oil standards.

Construction

The stainless steel construction ensures an exceptional stable bath temperature which further is improved by an ingenious stirring mechanism with baffle plates. All wetted parts are made of stainless steel. This provides resistance against all usual bath fluids. The bath is fitted with four wheels, two of them having brakes. The cover of the bath is on request and has to be ordered separately. There's a wide range of possibilities such as round ø51 mm openings, ø61 mm openings, the use of a leveling platform or other square or round openings with other dimensions. A bath overflow outlet protects against overflow when bath fluid expands or raises due to submersion of parts.

Agitation

A vane type stirrer with maintenance free bearings moves the bath fluid past a special heater ensuring optimal bath temperature control and ensuring excellent uniformity inside the TV16000.

Span lies from 10° C up to $+60^{\circ}$ C (50° F / 140° F).

Safety

The bath conforms to CE regulation. Further the bath is equipped with a mechanical over temperature device which trips when in case of malfunction the bath exceeds the preset maximum temperature. This feature guarantees safe around the clock operation.

Optional

The system overall accuracy is within ± 0.005°C based on measurement with water as bath fluid. This value is a min/max recorded over one hour. Standard deviation is better than ± 0.002°C under these conditions.

amson Instruments **Specification**

Master Viscometers and Viscosity Oil Standards

ASTM D2162

Fine adjustment and offset

After the temperature control is stable, the offset can be adjusted in the range of -5.00°C to +5.00°C, if necessary. The temperature can be set in 0.001°C resolution. A special offset offers ± 0.005°C resulting in setting of the set point temperature in ± 0.005°C.

Ambient condition

For proper cooling performance ambient temperature needs to be within the range of 18 .. 23°C.

Heating / cooling rate

Heating rate 12°C/hour Cooling rate 2.2°C/hour


Accuracy of control


Stability min / max over one hour measured in water, please see graphs on the next page @10 °C $min / max \pm 0.005$ °C

Standard deviation ± 0.0019 °C @20 °C $min / max \pm 0.003$ °C Standard deviation ± 0.0013 °C @40 °C

 $min / max \pm 0.003$ °C Standard deviation ± 0.0013 °C

Homogeneity ± 0.003 °C

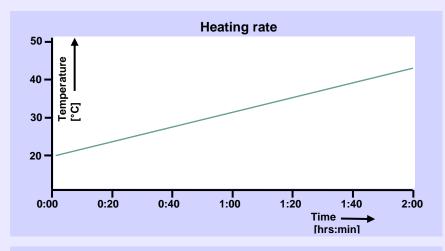

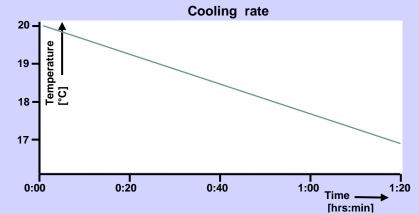






Tamson Instruments Specification sheet

Master Viscometers and Viscosity Oil Standards

ASTM D2162


@40 °CMin / max ± 0.003 °C
Standard deviation ± 0.0013 °C

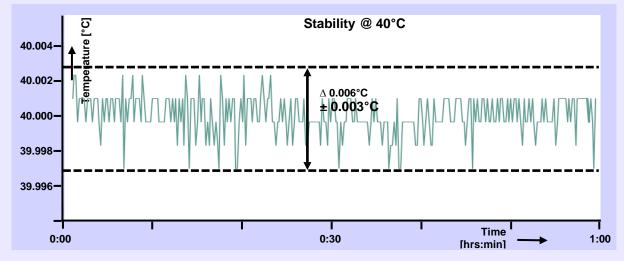

@20 °C

Min / max ± 0.003 °C


Standard deviation ± 0.0013 °C

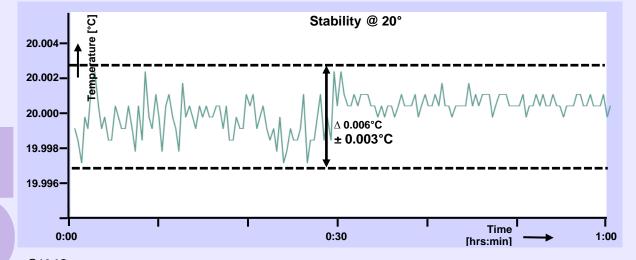

@10 °C

Min / max ± 0.005 °C

Standard deviation ± 0.0019 °C

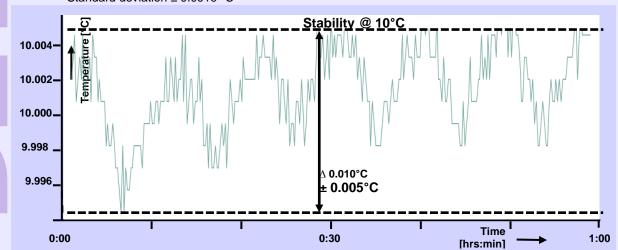

Tamson Instruments bv 🏻 🏗 31 105 22 43 73 🚖 sales@tamson.nl 🎓 www.tamson.com

Tamson Instruments Specification shee

Master Viscometers and Viscosity Oil Standards

ASTM D2162

P/N	Picture	Description
25T0594D		Master Ubbelohde viscometer with ISO 17025
25T0581P		calibration certificate size 0
25T0582P		Master Ubbelohde viscometer with ISO 17025
2510562F	₩	calibration certificate size 0B
25T0583P		Master Ubbelohde viscometer with ISO 17025
23103031		calibration certificate size 0C
25T0584P		Master Ubbelohde viscometer with ISO 17025
20100041		calibration certificate size 1
25T0585P		Master Ubbelohde viscometer with ISO 17025
2010000		calibration certificate size 1B
25T0586P		Master Ubbelohde viscometer with ISO 17025
		calibration certificate size 1C
25T0587P		Master Ubbelohde viscometer with ISO 17025
		calibration certificate size 2
25T0588P		Master Ubbelohde viscometer with ISO 17025
		calibration certificate size 2B
25T0589P		Master Ubbelohde viscometer with ISO 17025
		calibration certificate size 2C Master Ubbelohde viscometer with ISO 17025
25T0590P		calibration certificate size 3
		Master Ubbelohde viscometer with ISO 17025
25T0591P		calibration certificate size 3B
		Master Ubbelohde viscometer with ISO 17025
25T0592P		calibration certificate size 3C
		Master Ubbelohde viscometer with ISO 17025
25T0593P		calibration certificate size 4
		Master Ubbelohde viscometer with ISO 17025
25T0594P		calibration certificate size 4B
OFTOFOED		Master Ubbelohde viscometer with ISO 17025
25T0595P		calibration certificate size 4C
25T0596P		Master Ubbelohde viscometer with ISO 17025
2510596P		calibration certificate size 5
10T6035		Viscometer holder Master Ubbelohde
10T6030		Viscometer holder Master Cannon Fenske
On request		Uncalibrated master viscometers either Cannon
J 344001		Fenske or Ubbelohde.

Tamson Instruments Specification she

Master Viscometers and Viscosity Oil Standards

ASTM D2162

TV16000 accessories – options				
P/N	Picture	Description		
On request	○ —•	Top lid closed		
On request	000	Top lid ø51 mm openings		
On request	000	Top lid ø61 mm openings		

Alternative set-up TV7000DC + TLC15-5 with range from +15°C to +230°C and a temperature stability of \pm 0.007°C. For further information about the TV7000DC, please see specification sheet "TV7000DC".





